



PRINCIPLE AND GENERAL DESCRIPTION OF PIG 4MID[®] KIT (Ref. 4VDX-18K2)

The Pig 4MID[®] Kit (4VDX-18K2) is a sandwich ELISA assay that aims to detect and quantify proAKAP4 in boar sperm samples (in ejaculate, chilled or frozen semen in extender or in isolated spermatozoa). The Pig 4MID[®] Kit is composed of a 96-well microplate (12 x 8-well strips), and all the reagents and buffers required to run the assay. The proAKAP4 is a sperm-specific protein that is marker of sperm quality and male fertility in boars and main mammals (Ruelle et al. 2019 ; Sergeant et al. 2019 ; Delehedde et al. 2019 ; Carracedo et al. 2020 ; Ruelle et al. 2020 ; Griffin et al. 2020 ; Bastan and Akcay, 2021). Spermatozoa without proAKAP4 / AKAP4 are abnormal, immotile and infertile (Miki et al. 2002 ; Fang et al. 2019 ; Delehedde et al. 2019). Oxidative stress decreases proAKAP4 concentrations in spermatozoa (Nixon et al. 2019 ; Delehedde et al. 2019 ; Sergeant et al. 2020). Each 4MID[®] assay requires two major steps as described below in Part A and Part B of the Pig 4MID[®] Kit procedure. The proAKAP4 protein should be first released from spermatozoa using the Pig Spermatozoa Lysis Buffer and according to the types and/or sperm concentrations of the samples (ejaculate, semen in extenders or isolated spermatozoa) as described in Part A of the Pig 4MID[®] procedure. Then in Part B of the procedure, each prepared semen sample is loaded onto a 96-well microplate where a first antibody coated onto the bottom of the microplate will capture the proAKAP4 protein contained in the samples and will then be identified using a second antibody that is covalently coupled to horseradish peroxidase (sandwich method). A Substrate Solution is then added to each well and different color levels (blue) will appear proportionally to the concentrations of the proAKAP4 present in each semen sample. The color reaction is then stopped by adding the Stop Solution and the color intensity (yellow) is measured by spectrophotometry (reading at 450 nm). A positive control is included in each Pig 4MID[®] Kit. Ready-to-use Standard Solutions are provided to run a Linear Standard Curve allowing the calculation of proAKAP4 concentrations in boar semen samples. To compare animals or ejaculates or doses, we always recommend to express proAKAP4 concentrations in ng per 10 million of spermatozoa (ng/ 10 M of spz). Therefore, the concentrations in number of spermatozoa of the samples should be known before running the assay or be performed independently of the Pig 4MID[®] assay.

Always read the following instructions carefully before use

TABLE OF CONTENTS

I.	REAGENTS AND MATERIALS (INCLUDED)	2
II.	REAGENT AND MATERIAL REQUIRED - NOT INCLUDED	2
III.	STORAGE INFORMATION	2
IV.	GENERAL INSTRUCTIONS OF USE	2
A.	Preparation of the Semen Samples	3
a)	<i>Ejaculate or sample with sperm concentration above 50 million of spermatozoa per mL</i>	3
b)	<i>Semen in extenders or any sample with sperm concentration comprised between 10 to 50 million of spermatozoa per mL</i>	3
c)	<i>Isolated spermatozoa</i>	3
B.	Dosage of proAKAP4 Concentrations Using the Pig 4MID[®] Kit	4
C.	Calculation of Results – ProAKAP4 Concentration Determination	6
V.	PRACTICAL ADVICE AND CAUTIONS	7
VI.	REFERENCES	7
VII.	RELATED PRODUCTS	8



I. REAGENTS AND MATERIALS (INCLUDED)

- R1 - Capture Antibody Coated Microplate of 96-well (12 x 8-well strips)
- R2 - 1 Bottle of 10x Washing Buffer Solution (30 mL)
- R3 - 1 Bottle of 1x Dilution Buffer (30 mL)
- R4 - 7 Vials of Ready-to-Use Standard Solution (0.7 mL per vial)
- R5 - 1 Bottle of 1x Pig Spermatozoa Lysis Buffer (30 mL)
- R6 - 1 Vial of Detection Antibody (0.1 mL)
- R7 - 1 Bottle of Substrate Solution (11 mL)
- R8 - 1 Bottle 1x Stop Solution (7 mL)
- R9 - 1 Vial of Positive Control (0.7 mL)
- R10 - 2 Adhesive Plate Sealers
- R11 - User Manual Instruction

II. REAGENT AND MATERIAL REQUIRED - NOT INCLUDED

- 96-well microplate reader measuring absorbance at 450 nm
- Aluminum foil
- Horizontal orbital microplate shaker (300 rpm)
- Multichannel pipette of 300 µL
- One glass bottle (for R2 dilution)
- One polypropylene tube of 15 mL (for R6 dilution)
- Pen
- Pipette tips
- Pipettes of 20 µL, 200 µL and 1000 µL
- Plastic reagent reservoirs
- Polypropylene microtubes of 1.5 mL (for sample preparation)
- Scissors
- Ultrapure or double deionized water
- Vortex

III. STORAGE INFORMATION

- The Pig 4MID[®] Kit should be stored at 4°C upon receipt.
- All reagents must be protected from intense light.

IV. GENERAL INSTRUCTIONS OF USE

- Before use, bring all reagents except the R6 vial **at room temperature (RT) for at least 30 minutes before running the assay**. The R6 Detection Antibody vial should always be kept at 4°C.
- Verify the absence of crystals in the R2 and the R5 bottles. In presence of crystals, gently agitate the solution until all crystals are completely dissolved. Then R2 and R5 solutions should be kept at ambient temperature.
- The R1 96-well Plate (12 x 8-well strips) is in a reusable aluminum foil pouch. The plate frame and unused strips can be placed back in the reusable foil pouch for later use.



A. Preparation of the Semen Samples

ProAKAP4 should be first extracted from spermatozoa flagellum by using a specific R5 Pig Spermatozoa Lysis Buffer.

Do not forget before any pipetting of semen sample (fresh raw ejaculate, frozen semen in straws or isolated spermatozoa) to **resuspend cells by gently shaking** the tube containing the semen as spermatozoa cells will pellet by sedimentation.

Please note that before lysing spermatozoa cells, the semen should always be brought to ambient temperature but **never kept on ice**.

a) *Ejaculate or sample with sperm concentration above 50 million of spermatozoa per mL*

1. In a 1.5 mL conic tube, add 190 μ L of R5 Pig Spermatozoa Lysis Buffer.
2. Resuspend spermatozoa cells in the semen by gently shaking the semen before pipetting for homogenizing the sample.
3. Add 10 μ L of semen to the R5 Pig Spermatozoa Lysis Buffer to reach a volume of 200 μ L.
4. Carefully vortex each tube during 1 min **continuously** (minimum speed of 3000 rpm). All tubes should be vortexed before step 5.

Remark: Improper vortexing will lead to negative or false quantification value.

5. Add 200 μ L of R3 Dilution Buffer each sample.
6. Vortex rapidly at maximum speed.
7. Keep the lysed samples at ambient temperature (17°C – 25°C), not more than 6 hours before use. Do not put on ice.

b) *Semen in extenders or any sample with sperm concentration comprised between 10 to 50 million of spermatozoa per mL*

1. In a 1.5 mL conic tube, add 160 μ L of R5 Pig Spermatozoa Lysis Buffer.
2. Resuspend spermatozoa cells in the semen by gently shaking the semen before pipetting for homogenizing the sample.
3. Add 40 μ L of semen to the R5 Pig Spermatozoa Lysis Buffer to reach a volume of 200 μ L in each tube.
4. Carefully vortex each tube during 1 min at maximum speed **continuously** (minimum of 3000 rpm). All tubes should be vortexed before step 5.

Remark: Improper vortexing will lead to negative or false quantification value.

5. Add 200 μ L of R3 Dilution Buffer in each sample.
6. Vortex rapidly at maximum speed.
7. Keep the lysed samples at ambient temperature (17°C – 25°C), not more than 6 hours before use. Do not put on ice.

c) *Isolated spermatozoa*

1. Resuspend the spermatozoa pellet (between 700000 and 350000 spermatozoa) with 150 μ L of the R5 Pig Semen Lysis Buffer to reach a final concentration of 3.5 million sperm per milliliter in the R5 Pig Semen Lysis Buffer, not at concentration lower than 2.5 million sperm per milliliter in the R5 Pig Semen Lysis Buffer.

For instance, add 200 μ L of the R5 Pig Semen Lysis Buffer to a pellet of boar semen containing 0.7 million of spermatozoa.



User Manual Instructions – R11

2. Carefully vortex each tube during 1 min at maximum speed **continuously** (3000 rpm or above). All tubes should be vortexed before step 3.
Remark: Improper vortexing will lead to negative or false quantification value
3. Add 200 µL of the R3 dilution buffer in each sample.
4. Vortex rapidly at maximum speed
5. Keep the lysed samples at ambient temperature (17°C – 25°C) not more than 6 hours before use. Do not put on ice.

B. Dosage of proAKAP4 Concentrations Using the Pig 4MID[®] Kit

1. Open carefully the reusable aluminum foil pouch containing the R1 Microplate using scissors.
2. Add 100 µL of each Standard Solution (vial n°R4-1 to vial n° R4-7) on the first 8-well strip **from A1 well to G1 well** of the R1 Microplate (see table below). Please note that the R4-7 Standard Solution is the **Negative Control**. Always use a new tip for each different solution.

Standard vial n°	R4-1	R4-2	R4-3	R4-4	R4-5	R4-6	R4-7
ng / mL of proAKAP4	150	75	37.5	18.75	9.4	4.7	0
Well position	A1	B1	C1	D1	E1	F1	G1

Always follow this plate scheme if using the calculation datasheet provided by 4BioDx[®] services.

3. Add 100 µL of the **R9 Positive Control** in H1 well of the first strip of the R1 Microplate.
4. Then add **150 µL of each semen sample as prepared in Part A**. Up to 88 samples can be analyzed on the R1 Microplate from **well A2 to well H12** (from strip 2 to strip 12).

Remark 1: Vortex rapidly each semen sample before loading

Remark 2: Always use a new tip for each different semen sample.

5. Cover the R1 Microplate with one R10 Adhesive Plate Sealer and incubate the R1 Microplate for 2 hours at room temperature on a horizontal shaker with gentle agitation (300 rpm)
6. During the incubation time:

- a. Prepare the Detection Antibody: Dilute the content of the R6 Detection Antibody tube in 11 mL of R3 Dilution Buffer in a 15 mL tube. Close the tube and mix the Detection Antibody Solution by reversing the 15 mL tube several times. Keep this solution at ambient temperature until Step 9.

Remark 1: Before dilution in R3, please carefully hit the bottom of the R6 tube vertically or rapidly centrifuge to pull-down any liquid in the cap of R6 tube.

Remark 2: After dilution in R3, the Detection Antibody Solution should be used within one month and kept at 4°C.

- b. Prepare the Washing Solution: Add 30 mL of the R2 into 270 mL of ultrapure water (or double deionized water) in a clean glass bottle. Gently agitate the solution, avoid foaming and keep at ambient temperature.



7. At the end of incubation time, remove the R10 Adhesive Plate Sealer and eliminate the samples by reversion of the R1 Microplate (or aspirate when using an automatic microplate washer).
8. Transfer some Washing Buffer in a plastic reagent reservoir. Wash each well by adding **300 µL of R2 Washing Buffer 1x Solution** using the multichannel micropipette. Then discard the Washing Solution by reversion of the R1 Microplate. Repeat two times more. Please tap down gently the R1 Microplate on an absorbent dry paper to remove residual R2 Washing Buffer droplets between each washing step and before loading the R6 Detection Antibody
Remark: Results of the assay will markedly be influenced by the proper performance of the washing steps.
9. Transfer all the Detection Antibody Solution in a plastic reagent reservoir. Add **100 µL of Detection Antibody Solution** prepared in step 6 to each well of the 96-well Microplate (R1) using the multichannel micropipette.
10. Cover the R1 Microplate with a new R10 Adhesive Plate Sealer and incubate the R1 Microplate **for one hour** at room temperature with gentle agitation (300 rpm).
11. Remove the R10 Adhesive Plate Sealer and eliminate the Antibody Solution by plate reversion (or aspiration when using an automatic microplate washer).
12. Transfer some Washing Buffer in a plastic reagent reservoir. Wash each well by adding **300 µL of R2 Washing Buffer 1x Solution** using the multichannel micropipette (or an automatic microplate washer). Then discard the Washing Solution. Repeat two times more. Please tap down gently the microplate on an absorbent dry paper to remove residual Washing Buffer droplet between each washing step and before adding the R7 Substrate.
Remark: Results of the assay will markedly be influenced by the proper performance of the washing steps.
13. Transfer the Substrate Solution in a plastic reagent reservoir Add **100 µL of R7 Substrate Solution** to each well using the multichannel pipette.
Remark 1: Keep carefully away from light with an aluminum sheet.
Remark 2: Please note that R7 Substrate Solution and R8 Stop Solution (step 15) should be added to the wells in the same order and with the same time interval.
14. Protect from light and incubate the R1 Microplate under gentle agitation (300 rpm) for **15 minutes** at room temperature.
Remark: The incubation must not exceed 30 minutes.
15. Transfer the Stop Solution in a plastic reagent reservoir. Add **50 µL of R8 Stop Solution** to each well using the multichannel pipette and place the R1 Microplate 2 minutes on an orbital shaker at 300 rpm to mix well before reading the R1 microplate.
16. Determine the optical density (OD) of the samples using a microplate reader set to 450 nm.
*Remark: Always perform the measure **immediately** after adding the Stop Solution.*



C. Calculation of Results – ProAKAP4 Concentration Determination


A Standard Curve must be performed for each analysis to calculate proAKAP4 concentrations in boar semen. You can either do the Standard Curve a) using the Calculation Data Sheet provided for each sperm preparation protocol (part A) or b) do it manually yourself using software such as Excel.

Remark: always take in account the dilutions you have made when considering sperm concentration of the semen samples.

a) Using the Calculation Data Sheets provide by 4BioDx[®] services (upon request at contact@4biodx.com). The Calculation Data Sheets have been designed to automatically draw the Standard Curve from your optical density values (OD) and calculate the proAKAP4 concentrations first in ng /mL, then in ng/10 millions of spermatozoa. First select the calculation data sheet according to the sample preparation procedure (a, b or c) you have used in Part. A. You will then need to fill the blue cases with OD obtained from the plate reader and to follow the indicated steps on the sheet to draw the curve and then the sheet will automatically calculate the concentrations of proAKAP4 in ng/mL. By adding the concentration of spermatozoa in millions per mL of the semen samples, or the number of spermatozoa in the case of isolated spermatozoa (blue column of the step 5), you will obtain the concentrations of proAKAP4 in ng / 10 millions of spermatozoa. You may then copy and paste the values, tables and graphs on your own report files. Please note that hitting the step 6 button will generate a new sheet with summary of the results.

b) Manually, you will need to first subtract the optical density obtained from value G1 to the optical density of each standard (A1 to G1) and each sample (A8 to H12) that you obtained from the plate reader. Then create a Standard Curve by reporting the data (optical density values) on an Excel Spreadsheet. Express the optical densities (OD) in abscissa in the function of the proAKAP4 quantities in ng/mL in ordinate and then generate a two-degree polynomial regression equation (ProAKAP4 [conc] = a x (OD)² + b x (OD)) using the data analysis tools of an Excel Spreadsheet.

ProAKAP4 in (ng / mL)	Optical density (OD) values
150	OD A1
75	OD B1
37.5	OD C1
18.75	OD D1
9.4	OD E1
4.7	OD F1
0	OD G1

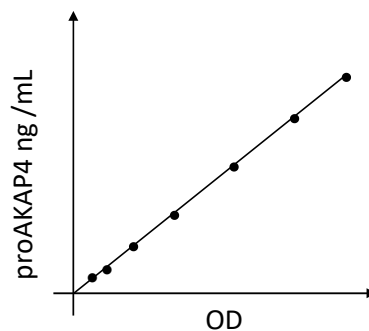

Then to **express the proAKAP4 concentrations in ng/mL**, calculate the concentrations using a two-degree polynomial regression equation.

To compare animals or ejaculates or doses, we recommend to express proAKAP4 concentrations **in ng per 10 million of spermatozoa (ng/10 M of spz)**.

To express results in ng of proAKAP4 per 10 million of spermatozoa, the formula will be:
 Conc. proAKAP4 ng/ 10 M = (conc. proAKAP4 in ng/mL / conc. spz M/mL) x 10 x dilution factor* x (2/3)
 (*use the dilution factor of 40 and 13.3 for the sample preparation procedure a) or b), respectively).



In the case of isolated spermatozoa: you should first calculate the concentration of spermatozoa in Million per mL (M/mL). To do so, multiply by 5 the number of spermatozoa in the pellet (between 700 000 and 500 000 of spermatozoa at step 2a of the semen sample preparation part).

Express then the results in ng of proAKAP4 per 10 million of spermatozoa (ng/ 10 of spz) with the following formula: Conc. proAKAP4 in ng/ 10 M of spz = (conc. proAKAP4 in ng/mL / conc. spz M/mL) x 10 x 0.66

V. PRACTICAL ADVICE AND CAUTIONS

- For research use only purposes and not for use for diagnostic purposes.
- Always follow good laboratory practices.
- Always follow the plate scheme indicated in part C when using the Calculation Datasheet provided by 4BioDx[®] services.
- After dilution in R3, the Detection Antibody Solution should be used within one month and kept at 4°C.
- The use of a multichannel pipette is mandatory to ensure the timely delivery of liquids.
- Use the supplied reagents as an integral unit prior to the expiration date.
- The R7 Substrate Solution can be irritating for the skin.
- The R8 Stop Solution can be harmful in case of ingestion and could lead to irritation when in contact with the skin.
- To avoid distortions due to differences in incubation times, R7 Substrate Solution and R8 Stop Solution should be added to wells in the same order and with the same time interval.
- Do not expose the R7 Substrate Solution to light or oxidative substances.
- Use only reagents from the same Pig 4MID[®] Kit.
- Any variation in ambient temperature, pipetting, washing method or incubation time can cause variation in optical density results.
- Always use new tips for each different solution and for each sample.
- Observe all federal, state, and local regulations in terms of waste disposal.

VI. REFERENCES

Bastan I and Akcay E (2021) Quality assessment of frozen bull semen with the precursor A-kinase anchor protein 4 biomarker. *Andrologia* e14164.

Xu Y, Han Q, Ma C, Wang Y, Zhang P, Li C, Cheng X and Xu H (2021) Comparative proteomics and phosphoproteomics analysis reveal the possible breed difference in Yorkshire and duroc boar spermatozoa. *Front Cell Dev Biol.* Vol. 9:652809. doi: 10.3389/fcell.2021.652809. eCollection 2021.

Blommaert D, Sergeant N, Delehedde M, Donnay I, Lejeune JP, Franck T and Serteyn D (2021) First results about ProAKAP4 concentration in stallion semen after cryopreservation in two different freezing media. *Cryobiology.* Vol. 102:133-135.

Carracedo S, Loyens A, Eddarkaoui S, Serteyn D, Malo C, Skidmore L, Briand-Amirat L, Barbotin AL, Mauraage CA, Delehedde M and Sergeant N (2020) The sperm specific proAKAP4 polypeptide exhibited conserved functions, localizations and metabolism among mammals. *Animal Reproduction Science.* Vol. 220 :106448

Riesco M, Anel-Lopez L, Neila-Montero M, Palacin-Martinez C, Montes-Garrido R, Alvarez M, de Paz P, Anel L (2020) ProAKAP4 as Novel Molecular Marker of Sperm Quality in Ram: An Integrative Study in Fresh, Cooled and Cryopreserved Sperm. *Biomolecules.* 10(7):1046.

Sergeant N, Carracedo S, Blommaert D, Aubry S, Jouy N, Lejeune JP, Franck T, Mauraage CA, Serteyn D, Buée L and Delehedde M (2020) Proteolysis of proAKAP4 in semen as a regulatory sensor of sperm quality and functionality. *Andrology.* Vol. 8 (Suppl. 1):44-45.



Ruelle I, Sergeant N, Bencharif D, Charreaux F, Thorin C, Michaud S, Dordas-Perpinyà M, Jouy N, Audry S, Maurage CA, Delehedde M and Briand-Amirat L (2020) ProAKAP4 concentrations in semen as a predictive tool of bull fertility: A preliminary study. *Reproduction, Fertility and Development*. 32(1):145.

Griffin R, Swegen A, Baker M, John Aitken R, A Skerrett-Byrne D, Silva Rodriguez A, E Martín-Cano F, Nixon B, J Peña F, Delehedde M, Sergeant N and Gibb Z (2020) Mass spectrometry reveals distinct proteomic profiles in high- and low-quality stallion spermatozoa. *Reproduction*. 160(5):695-707.

Nixon B, Bernstein I, Cafe SL, Delehedde M, Sergeant S, Eamens AL, Lord T, Dun MD, De Iuliis GN and Bromfield EG (2019b) A Kinase Anchor Protein 4 is vulnerable to oxidative adduction in male germ cells. *Frontiers in Cell and Developmental Biology* – Vol. 7:319.

Fang X, Huang LL, Xu J, Ma CQ, Chen ZH, Zhang Z, Liao CH, Zheng SX, Huang P, Xu WM, Li N, Sun L (2019) Proteomics and single-cell RNA analysis of Akap4-knockout mice model confirm indispensable role of Akap4 in spermatogenesis. *Developmental Biology*. 454(2):118-127.

Sergeant N, Briand-Amirat L, Bencharif D and Delehedde M (2019) The sperm specific protein proAKAP4 as an innovative marker to evaluate sperm quality and fertility. *Journal of Dairy & Veterinary Sciences*. Vol. 11:01-19.

Ruelle I, Charreaux F, Bencharif D, Thorin C, Michaud S, Schmitt E, Sergeant N, Briand-Amirat L and Delehedde M (2019) Assessment of the sperm specific protein proAKAP4 as a marker to evaluate sperm quality and fertility in Holstein bulls. *Rev. Bras. Reprod. Anim.* Vol.43(2):472.

Delehedde M, Carracedo S, Selleslagh M, Eddarkaoui S, Amirat-Briand L and Sergeant N (2019) ProAKAP4 polypeptide as a biomarker of sperm functionality and male fertility disorders. *Int J Gynecol and Reprod Sci*. Vol. 2(1):13-19.

Delehedde M, Demouveaux B, Remy G, Selleslagh M, Dewulf Q, Desseyn JL, Moreau M, Gosset P, Muriel Pichavant M and Sergeant N (2019) ProAKAP4 concentrations as an indicator of good spermatogenesis and sperm quality under oxidative stress conditions. *Andrology*. Vol. 7(1):86.

Peddinti D, Nanduri B, Kaya A, Feugang JM, Burgess SC and Memili E (2008) Comprehensive proteomic analysis of bovine spermatozoa of varying fertility rates and identification of biomarkers associated with fertility. *BMC Systems Biology*. Vol. (2)19:1-13.

Miki K, Goulding EH, Fulcher KD, Eddy E, Willis WD and Brown PR (2002) Targeted disruption of the AKAP4 gene causes defects in sperm flagellum and motility. *Developmental Biology*. Vol. 248:331-342.

VII. RELATED PRODUCTS

Reference	Designation	Specificity
4VDX-18K2	Pig 4MID® Kit	Pig proAKAP4
4VDX-18K3	Horse 4MID® Kit	Horse / Donkey proAKAP4
4VDX-18K4	Bull 4MID® Kit	Bull proAKAP4
4VDX-18K5	Dog 4MID® Kit	Dog proAKAP4
4VDX-18K6	Rabbit 4MID® Kit	Rabbit proAKAP4
4VDX-18K7	Ram 4MID® Kit	Ram proAKAP4
4BDX-18K8	Mouse 4MID® Kit	Mouse / Rat proAKAP4
4VDX-18K9	Goat 4MID® Kit	Caprine proAKAP4
4VDX-19K10	Cat 4MID® Kit	Cat proAKAP4 (feline)
4VDX-19K11	Camel 4MID® Kit	Camelids proAKAP4



Distribuito in ITALIA da
Li StarFish S.r.l.
 Via Cavour, 35
 20063 Cernusco S/N (MI)
 telefono 02-92150794
 info@listarfish.it
 www.listarfish.it

FOR RESEARCH USE ONLY PRODUCT
NOT FOR USE IN DIAGNOSTIC PROCEDURE